

SAFETY FEATURES

To prevent entrapment the buddy-system must be used.

Two individuals must be present each time the bulk sterilizer is started.

The inside of the autoclave is observed to confirm that no one is inside prior to activating.

Both individuals will remain outside the autoclave to ensure that no one is inside as the door is closed and the sterilizer cycle is being activated.

By design the door to the sterilizer only closes very slowly to reduce the possibility of entrapment.

Emergency Shut-Off (red button) located on both sides of the autoclave.

Emergency Electrical Cut-Off located to left of unit on prep-side.

SC #259
Eff 7/19